CENTRA

CENTRALIZED WATER




CENTRA-S/R 200

The CENTRA has revolutionised the way that large volumes of pure water are produced, stored and distributed. Instead of a central lab system, comprising of an untidy grouping of components and pipework, you can now have just one integrated system box. The CENTRA-S/R 200 are complete water purification, storage, control and distribution systems featuring a 200 liter per hour reverse osmosis module and 0.2 µm filter.

- The compact design provides more flexible installation options for new buildings and refurbishments. The CENTRA's small footprint means that it can be placed closer to the laboratory, avoiding the negative cost and design implications of long pipe work loops
- Reliable, continual supply of pure water using unique access controls, leak detection systems and full AV alarms with optional building management system (BMS) connectivity
- Optimized inorganic water quality (RDS) through use of in-line purification technologies. Recirculated water is UV treated, filtered and (if fitted) can be improved through deionization
- 200 l/hr of purified water available at up to 30 l/min from a distribution loop. A wide range of water purities are possible from RO permeate up to 18.2 M Ω -cm Type I purity
- Low microbial counts achieved through vent filtration, spray-ball water distribution and smooth internal surface in the reservoir combined with UV oxidation and 0.2 µm filtration in the loop

Process Flow CENTRA S/R 200

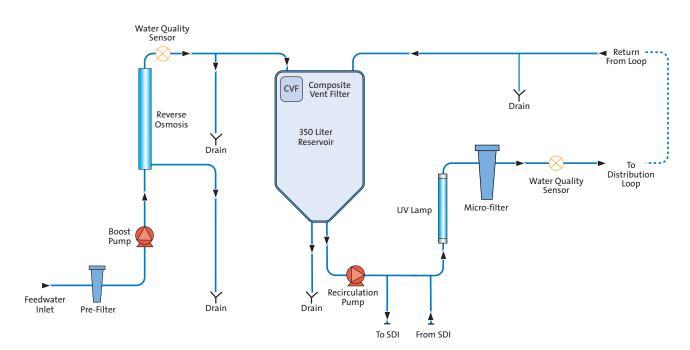





CENTRA

CENTRALIZED WATER

Treated Water Specifications

Model	Voltage/freq.	S/R 200	S/R 200 HFV	S/R 200 HFR
Flowrate (exit from unit) - I/min	230/50	16 (4.2)	30 (8)	30 (8)
(USG/min)	115/60	18 (4.8)	38 (10)	38 (10)
Daily usage - typical		Up to 5000 liters (1321 USG)	Up to 5000 liters (1321 USG)	Up to 5000 liters (1321 USG)
Daily usage - maximum		Up to 6000 liters (1585 USG)	Up to 6000 liters (1585 USG)	Up to 6000 liters (1585 USG)
Model	S200		R200	
Inorganics	RO Permeate ¹		Up to 18.2MΩ-cm²	
TOC - typical	<100 ppb1		<10 ppb³	
Bacteria - typical	<50 CFU/ml³		<5 CFU/ml³	
Particles	0.2 μm filtration		0.2 µm filtration	

¹ Feedwater dependent. ² Ion-exchange cylinder installed (Nuclear or Hypex grade). ³ System to be installed following ELGA LabWater installation guidelines and regularly sanitized.

Dimensions and weights

Dimensions	Height 1820mm (71.7in), Width 730mm (28.75in), Depth 890mm (35in)		
Shipping weight	180kg (396lb)	178kg (392lb)	
Operational weight	530kg (1168lb)	527kg (1162lb)	

Feedwater Requirements

Source Potable tap water as detailed below.

Note: It is essential that the feedwater be suitably conditioned. Please refer to the contaminants listed below and ensure that suitable pretreatment is included in the installation.

Contaminant	Measure	Range	Pretreatment	
Calcium	Ca ppm as CaCO ₃	<250	None ⁴	
		>250	Softener or use very low RO recovery ⁴	
Total chlorine	Cl ppm	<0.1	None	
		0.1 to 0.5	20 inch carbon block	
		>0.5	Cylinder of carbon sized correctly to obtain < 0.1ppm	
Silica	SiO₂ ppm	<30	None	
		>30	20 inch cartridge depth filter	
Fouling Index	FI	<10	None	
		10 - 20	20 inch cartridge depth filter	
		>20	Back washable media filter with a minimum flow rate of 20 l/mir	
Iron/manganese	Fe/Mn ppm	<0.05	None	
		>0.05	20 inch cartridge depth filter ⁶	
		>0.1	Back-washable pre-filter ⁶	
Organics	TOC ppm C	<2	None	
		2 - 3	20 inch carbon block ⁵	
		>3	Cylinder of carbon sized correctly for TOC demand ⁵	
Temperature	1 - 40°C (Recommended 15 - 25°C)			
Flowrate (maximum requirement @15°C)	20 l/min (5.3 GPM)			
Drain requirements (gravity fall with air gap)	45 l/min (12 GPM)			
Feedwater Pressure	4 bar (60 psi) maximum, 2 bar (30 psi) minimum.			

⁴ Check LSI, increase frequency of acid cleaning. ⁵ Increase frequency of alkaline cleaning. ⁶ Increase frequency of acid cleaning.

Electrical Requirements

Mains input	230V ac, 50Hz, 115V ac, 60Hz
System control voltage	24V dc
Power consumption during recirculation	2000VA
Electrical protection rating	20 amps
Reservoir level connection for control of additional ELGA products	Jack Plug 3.5mm DIN plug 6 way
Noise level during recirculation	<70dBA

ELGA LabWater

Tel: +44 (0) 1494 887500 Fax: +44 (0) 1494 887505

Email: info@elgalabwater.com Website: www.elgalabwater.com

ELGA® is the global laboratory water brand name of Veolia Water. VWS (UK) Ltd. Registered in England & Wales No. 327847 ©Copyright 2008 ELGA LabWater/VWS (UK) Ltd. All rights reserved. As part of our policy of continual improvement we reserve the right to alter the specifications given in this datasheet.